

Changing Gas Cylinders

Preparation

- For toxic gasses, take proper precautions
- Obtain correct regulator. Check CGA# (See SOP).
- Check special gas (SG) labelling on hood ports
- Check status of valves and connections

Operation

- Cylinders should be double strapped to wall while in use
- Never force a connection! Do not overtighten nut!**
- Delivery pressure: 5-10 psi
 - Gauge reading won't decrease in a closed system!**
- Leak check connections from cylinder to bubbler before use
- Never pressurize a Schlenk line — keep bubbler open