

Glovebox Failure Checklist

1. Gloves are pulled inside (*Low Pressure*)

- Turn off catalyst circulator
- Replace N₂ tank if empty
- Purge for 1 hour
- Turn on catalyst circulator

2. Gloves are very inflated (*High Pressure*)

- Open small antechamber and turn on vacuum
- Close small antechamber once pressure has stabilized

3. O₂ levels are >100 ppm

- Turn off catalyst circulator
- Purge for 1 hour
- During purge, check for leaks in antechamber and gloves
- Turn on catalyst circulator